

Tera-SED

Large area THz emitter

- Large-area terahertz emitter
- Low external bias voltages
- No external cooling required
- Superior interdigitated design

Overview


Efficient sources of Terahertz (THz) radiation are of great importance for a large variety of scientific and technological applications. Important key factors are a large bandwidth and a high THz electric field amplitude. This usually implies a large emitter area and in conventional concepts requires high bias voltages (several hundreds of volts) in order to achieve the necessary bias field strengths, typically in the kV/cm range. The key advantage of the Tera-SED is the unique combination of large active area and low bias voltage requirement.

Technical information


The **Tera-SED** is a planar large-area GaAs-based photoconductive emitter for impulsive generation of broadband THz radiation with a superior design. It features a novel interdigitated electrode MSM structure which allows for a large active area with kV/cm-bias fields between individual electrodes. Only low external bias voltages are required, eliminating the need for a pulsed, high-voltage supply. Its flexible scalable device technology makes **Tera-SED** an efficient and versatile THz emitter.

Tera-SED requires no external cooling. It comes attached to a metal holder fitting into 1-inch optics mounts. Ease of use and hassle-free alignment are key assets.

Tera-SED is available in two different versions: 10x10 mm² usable area intended for use with amplified fs-laser systems (pulse energy up to 300 μJ, saturating from ~10 μJ) and 3x3 mm² usable area best suited for use with fs-laser oscillators.

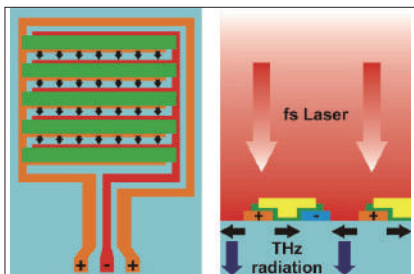

Tera-SED functional principle: with applied bias voltage the electric field direction is reversed between successive electrode fingers. An optically opaque metal layer is applied between every other finger pair such that optical excitation is only possible in areas exhibiting the same electric field direction. THz radiation emitted from the device thus interferes constructively in the far-field.

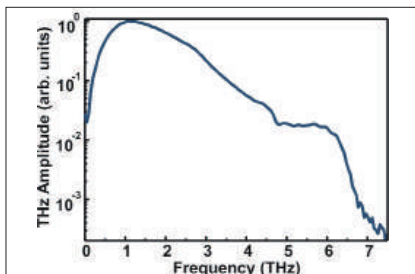

Tera-SED3 emission spectrum at 10 V DC bias and 0.7 nJ pulse energy, acquired with Laser Quantum GmbH (formerly Gigaoptics) ' High-Speed ASOPS technique.

脉动科技有限公司
中国代理商

总部地址：海淀区中关村东路84-8号 传真：010-62565117-803
 二层1-104单元，100190 邮箱：info@pulsepower.cn
 电话：010-62565117 官网：www.pulsepower.cn

上海办事处 021-60720971 深圳办事处 0755-26922729 成都办事处 17710735621



Key device characteristics

| | Tera-SED3 | Tera-SED10 |
|--|---------------------|-----------------------|
| Size | 3x3 mm ² | 10x10 mm ² |
| Peak emission frequency | 1.0-1.5 THz | |
| Spectral width (@-10 dB) | ~2.5 THz | |
| Pulse energy | 7.5 nJ | 10 μJ |
| Max.opt. excitation intensity | 8 W/mm ² | |
| Max.opt. excitation power | 650 mW | |
| Optical excitation wavelength | 700-850 nm | |
| Pulsed THz field amplitude: | < 5kV/cm | |
| Low bias voltage (DC or switchable) ¹ | 1-30 V | |
| Bias modulation frequency | DC to 100 kHz | |
| Duty cycle ² | 5% to 100% (cw) | |

¹ max. bias voltage depends on active area and type of usage.

² max. duty cycle depends on active area and type of usage.

Electrical performance parameters*

| | Vbias | duty cycle | THz field amplitude |
|---------------------|------------------|------------|---------------------|
| Tera-SED3 | up to 10 V | CW | 100 V/cm (@10 V) |
| 7.5 nJ pulse energy | 10 V-20 V | 50% | 200 V/cm (@ 20 V) |
| Spotsize | 300 μm max. 30 V | 10% | 300 V/cm |
| Tera-SED10 | up to 5 V | CW | 1000 V/cm (@5V) |
| 10 μJ pulse energy | 5 V-20 V | 50% | 2000 V/cm (@20 V) |
| Spotsize 2 to 6 mm | max. 25 V | 5% | 5000 V/cm |

* Typical values only. Actual field strengths may vary depending on specific experimental conditions.



LASER QUANTUM LTD

tel: +44 (0) 161 975 5300
email: info@laserquantum.com
web: www.laserquantum.com

LASER QUANTUM INC

tel: +1 510 210 3034
email: info@laserquantum.com
web: www.laserquantum.com

LASER QUANTUM GmbH

tel: +49 7531 368371
email: info@laserquantum.com
web: www.laserquantum.com